

# Shock sensor alarm

## Software installation guide

### **Contact Information**

Email: [info@laboratories.cz](mailto:info@laboratories.cz)

Web: [laboratories.cz](http://laboratories.cz)



## Contents

Driver installation.....	3
Shock Alarm software installation.....	6
Shock Alarm Supervisor software installation.....	9



## Driver installation

1. Connect sensor device to USB port on your computer and wait. "No driver found" message appears with "Shock Sensor Laboratories" name.




2. Open the "Device Manager" window. In "Other devices" section, click on "Shock Sensor Laboratories" and choose "Update Driver Software...".






3. Click on "Browse my computer for driver software".

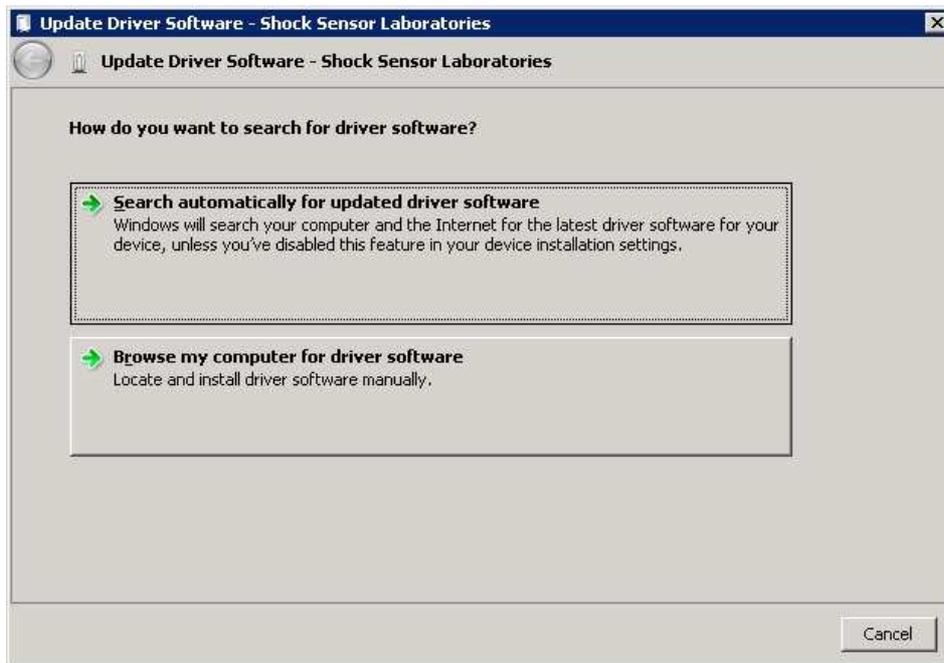

4. "Browse" to the folder where driver files are stored.

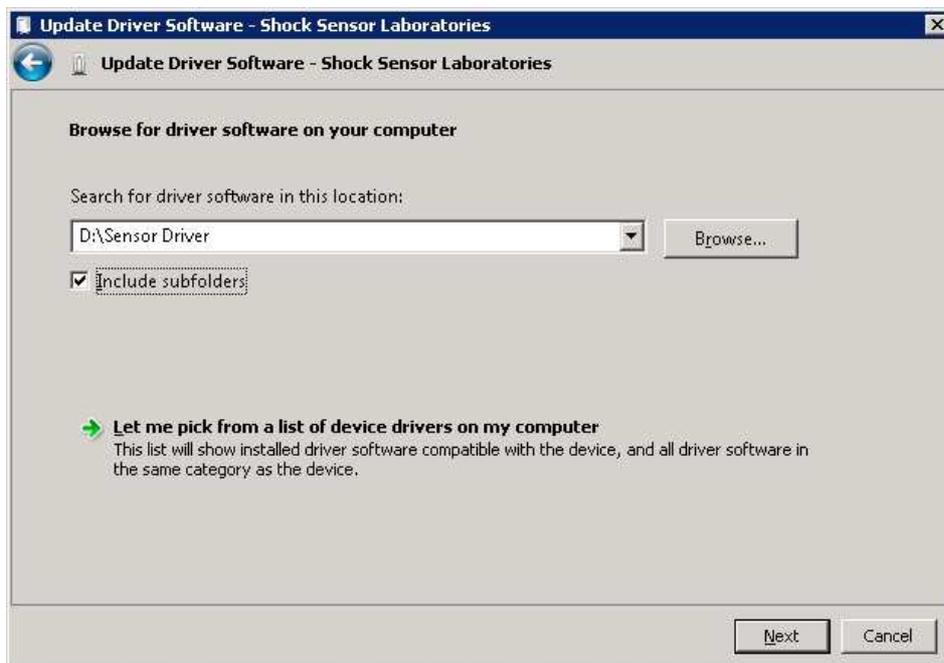

5. Click "Next".



6. Click "Close".



7. The Sensor should appear in "Universal Serial Bus controllers" section in the "Device manager".



8. Device driver was successfully installed.




## Shock Alarm software installation

**Prerequisites:** .NET Framework 2.0 installed on the target computer.


1. Run the ShockAlarm installation file.




2. Welcome window appears. Click "Next".

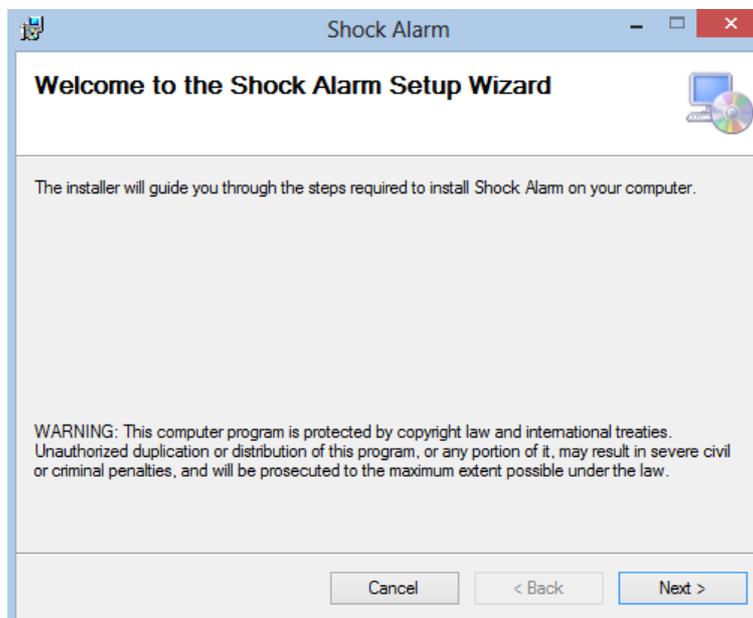


3. In the "Select Installation Folder" window set an installation folder and click "Next".

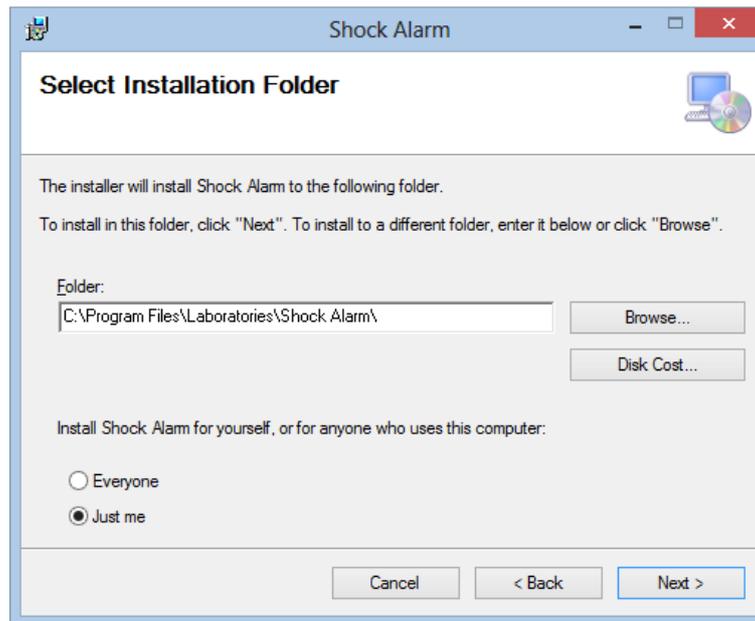


4. In the "Confirm Installation" window click "Next".

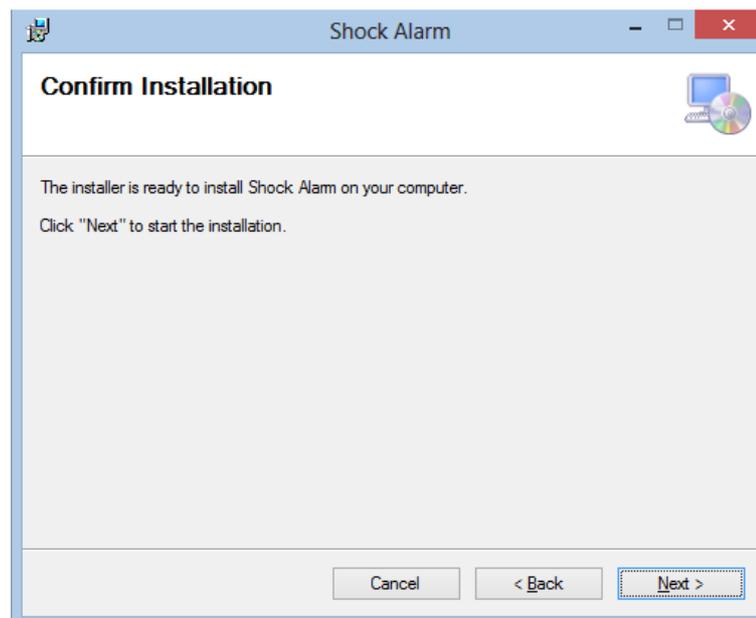

5. If the "User Account Control" message appears, click "Yes".

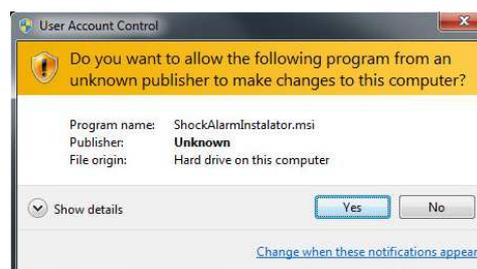

6. The "Installation Complete" window appears. Click "Close".

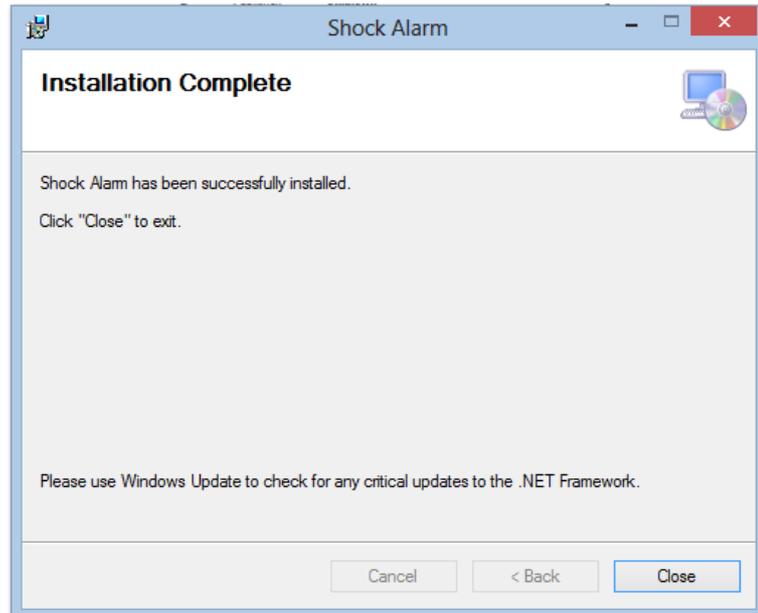

7. Run the Shock Alarm application from Windows desktop.






## Shock Alarm Supervisor software installation

**Prerequisites:** .NET Framework 4.0 installed on the target computer.


1. Run the ShockAlarm Supervisor installation file.

ShockAlarmSupervisorInstalator.msi

2. Welcome window appears. Click "Next".

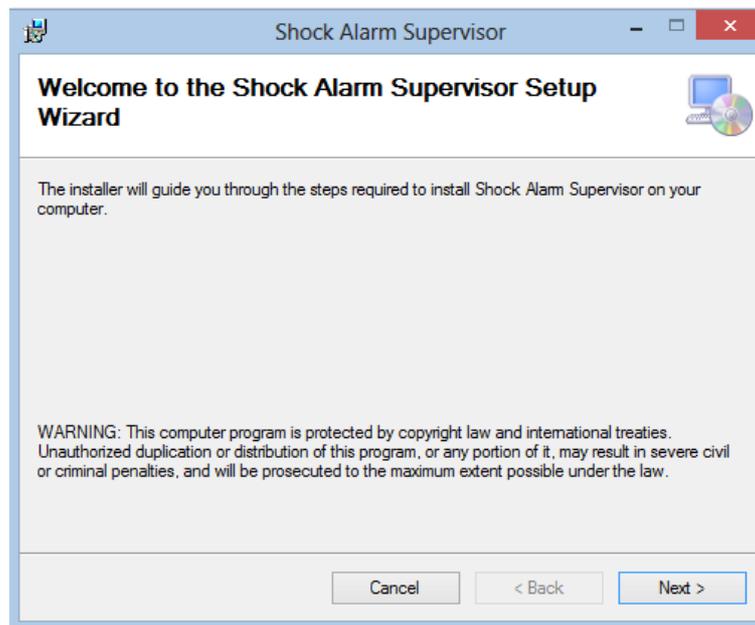


3. In the "Select Installation Folder" window set an installation folder and click "Next".

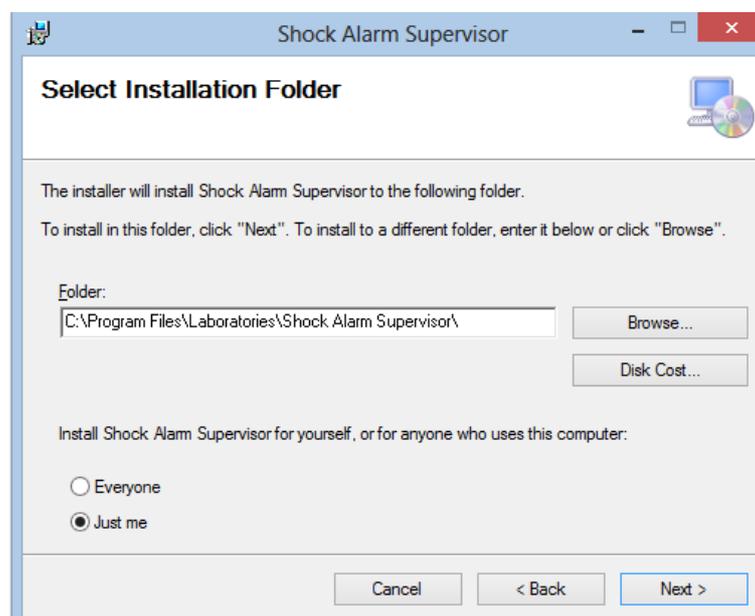





4. In the “Confirm Installation” window click “Next”.

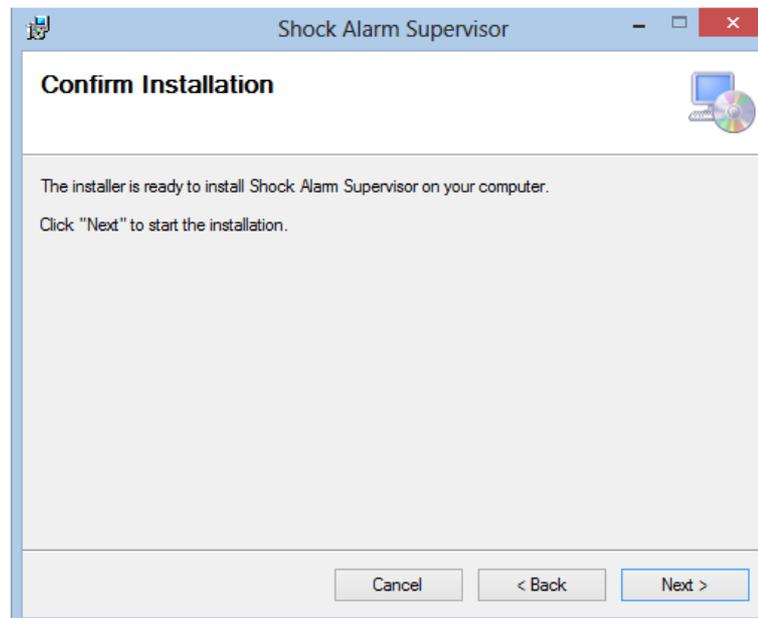

5. If the “User Account Control” message appears, click “Yes”.

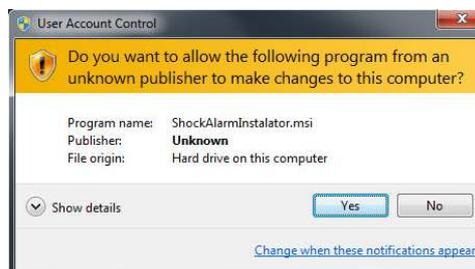

6. The “Installation Complete” window appears. Click “Close”.

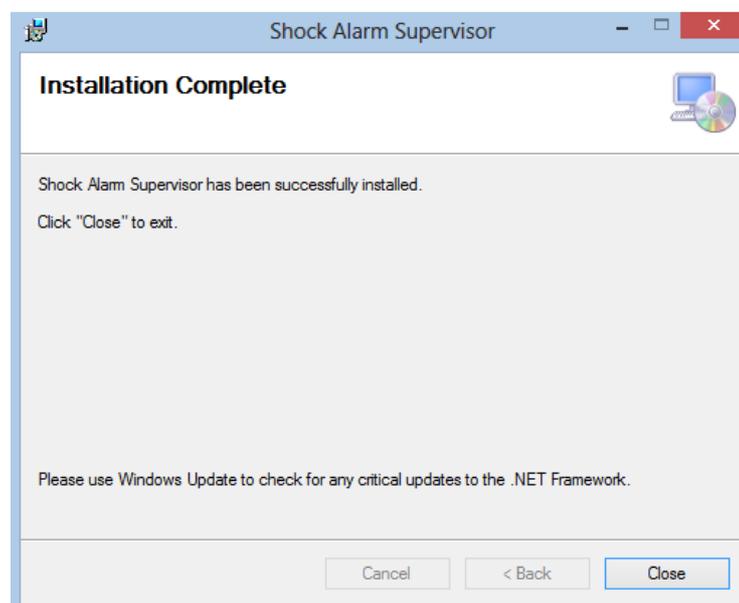



7. Run the Shock Alarm Supervisor application from Windows desktop.



**In case of any problems, please contact technical support Laboratories s.r.o.  
E-mail: [info@laboratories.cz](mailto:info@laboratories.cz)**